

**TO REPLACE PILOT LIGHT**

**#101409 IN EA-1**

1. Remove dashpot and piston
2. Remove counter weights
3. Remove weight rack and weights
4. Remove pan and tape doors closed
5. Gently tip balance back on its back
6. Remove bottom cover
7. Remove wire ties and disconnect pilot light
8. Push light out of console
9. Push new light into console and reconnect, replace wire ties
10. Retrace steps 6, 5, 4, 3, 2 & 1

**\*NOTE: Make sure wires do not interfere with counter.**